

SquidTM

Full Range Pipette (1 - 1000 μ L)

Quick Start Guide

2-APR-2026

DeNovix[®]

Active Pipetting Mode

See User Guide for pipetting mode instructions.


Tip Status

Action Wheel

Hold for 0.5 seconds to wake from Sleep Mode.

Press to aspirate, dispense or select.

Turn to scroll or adjust volume.

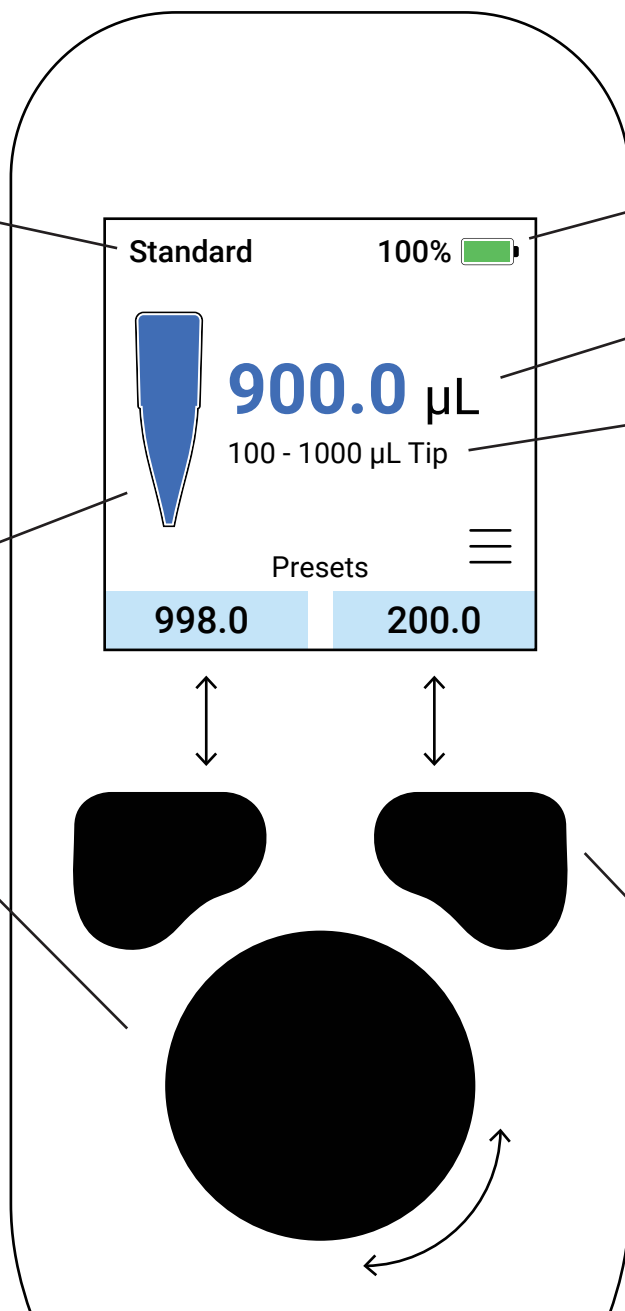

Battery

Volume

Tip Mode

Side Button

Access menu or return to previous screen.

Press and hold to put into Sleep Mode.

Left and Right Buttons

Press to activate Presets.

Hold to edit value.

Squid™ Full Range Pipette

Squid™ is a 1 - 1000 μL electronic pipette that uses two tip sizes for different volume ranges.

- Large Tip Mode: 100 - 1000 μL
- Small Tip Mode: 1 - 200 μL


What's in the Box?

- Squid™ Full Range Pipette (1 - 1000 μL)
- Charging Cable and Charger
- Certification
- O-Ring Grease
- Extra O-Rings

Getting Started

1. Remove Squid Pipette from packaging.
2. Hold Action Wheel for 0.5 seconds to wake from Sleep Mode.
3. Use the Side Button to select a pipetting mode or adjust settings.

For detailed instructions, please refer to the User Guide (QR code on back).

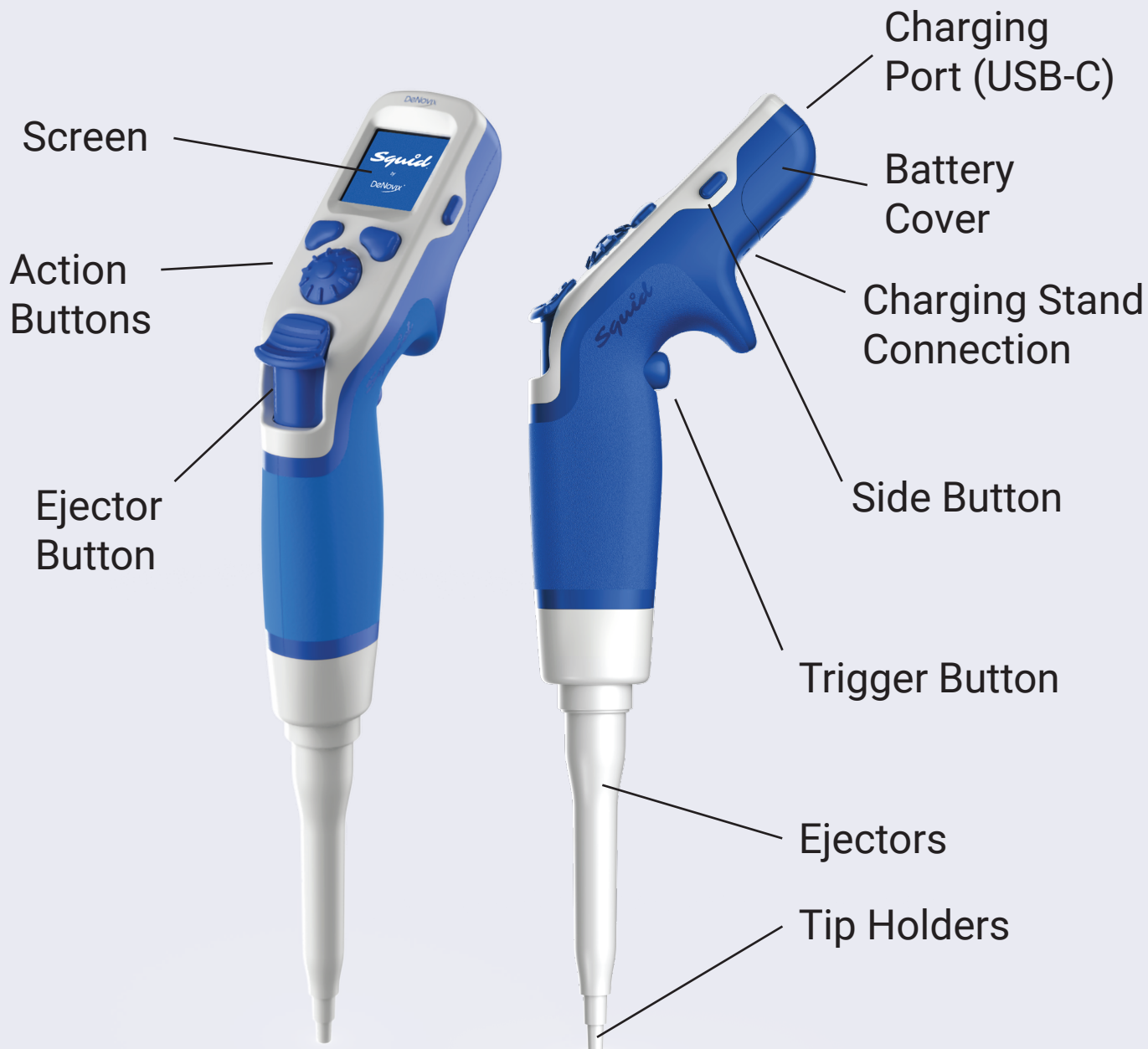

Using Squid Tip Modes

Large Tip Mode

100 - 1000 μ L



Press Ejector



Large Tip Ready

Small Tip Mode

1 - 200 μ L



Press Trigger Button



Small Tip Ready

Important: After ejecting, press Trigger Button to return to Small Tip Mode.



User Guide



Request Support
techsupport@denovix.com



Give Us Feedback



#SquidSquad