

# User Guide

## Orangu™

### Cell Counting Solution

Cat OR01

Protocol Version 2.0





## Contents

Product components .....	4
Storage .....	4
Equipment and materials required but not supplied with these reagents .....	4
Introduction and Assay Principle .....	5
Procedure .....	6
General protocol .....	6
Cell Proliferation and Cytotoxicity Assay .....	6
Notes on use of Orangu™ .....	8
Troubleshooting .....	9
References.....	9
Purchaser Notification .....	11

# Orangu™ Cell Counting Solution

## Product components

Table 1. Orangu™ range of products.

Product	Catalogue Number	Number of tests
Orangu™ Solution, 1 ml x 1	OR01-100	100
Orangu™ Solution, 5 ml x 1	OR01-500	500
Orangu™ Solution, 5 ml x 2	OR01-1000	1000
Orangu™ Solution, 5 ml x 6	OR01-3000	3000

## Storage

Store at 4°C. Protect from light. Orangu™ has a shelf life of at least six months from receipt.

## Equipment and materials required but not supplied with these reagents

- Absorbance microplate reader fitted with 450 nm filter
- 96-well plate (or multiwell plate of choice)
- CO<sub>2</sub> incubator
- Pipettes

## Introduction and Assay Principle

Orangu™ is a cell proliferation and cytotoxicity assay which utilizes WST-8, a water-soluble tetrazolium salt, which produces a formazan orange dye upon reduction in the presence of an electron mediator (Figures 1 and 2).

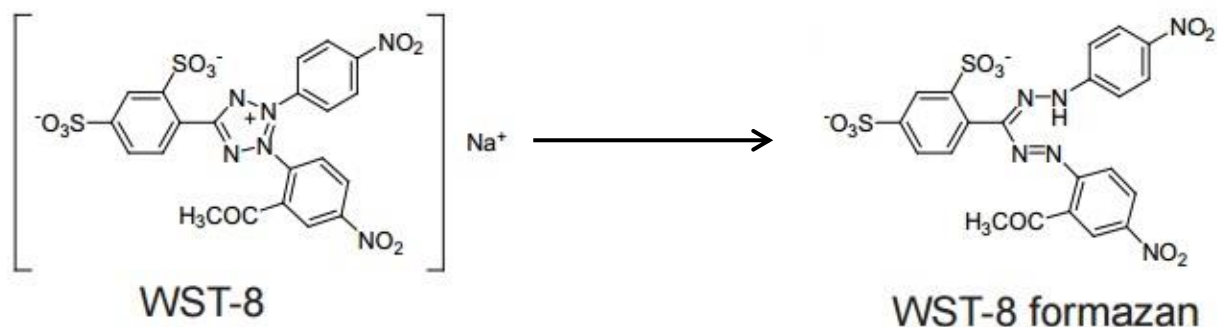


Figure 1. Structures of WST-8 and WST-8 formazan.

Orangu™ is a one-bottle solution; no premixing of components is required. Orangu™ provides a sensitive colorimetric assay for the determination of the number of viable cells in cell proliferation and cytotoxicity assays.

The amount of the formazan dye generated by dehydrogenases in cells is directly proportional to the number of live cells and the length of incubation.

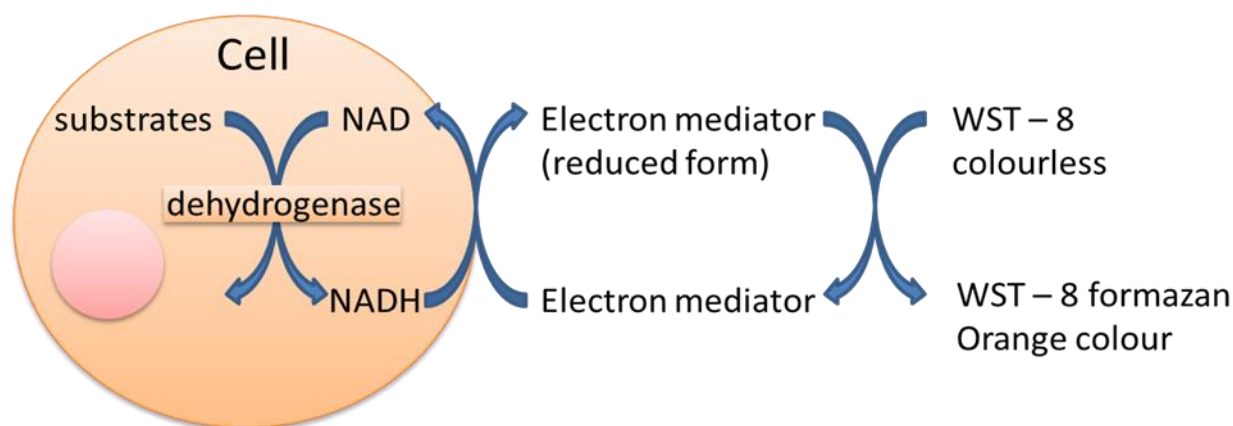


Figure 2. Principle of the cell viability detection with Orangu™.

Orangu™ data correlates very well with the [<sup>3</sup>H]-thymidine incorporation assay. Detection sensitivity using Orangu™ is higher than assays using other tetrazolium salts (such as MTT, XTT, MTS or WST-1).

## Procedure

If further culture of the cells is required following Orangu™ assay, please filter the solution through a 0.2 µm membrane to ensure sterility.

### General protocol

1. Inoculate cell suspension (100 µl/well) in a 96-well plate. Pre-incubate the plate in a humidified incubator (e.g. at 37°C, 5% CO<sub>2</sub>) to equilibrate.
2. Add 10 µl of Orangu™ solution to each well of the plate.  
**Do not introduce bubbles to the wells - they interfere with the absorbance reading.**
3. Incubate the plate for 1 - 4 hours in the humidified incubator.  
**More than 1 hour in the incubator allow a higher signal.**
4. Measure the absorbance at 450 nm using a microplate reader. To measure the absorbance later (within 24 hours), add 10 µl of 1% w/v SDS or 0.1 M HCl to each well, cover the plate and store it in the dark at room temperature.

### Cell Proliferation and Cytotoxicity Assay

1. Dispense 100 µl of cell suspension (5×10<sup>3</sup> cells/well) in a 96-well plate. Pre-incubate the plate for 24 hours in a humidified incubator (e.g. at 37°C, 5% CO<sub>2</sub>).
2. Add 10 µl of various concentrations of substances to be tested to the plate.
3. Incubate the plate for an appropriate length of time (e.g. 6, 12, 24 or 48 hours) in the incubator.
4. Add 10 µl of Orangu™ solution to each well of the plate. Be careful not to introduce bubbles to the wells, since they interfere with the absorbance reading.
5. Incubate the plate for 1 - 4 hours in the humidified incubator.  
**Incubation periods will depend on cell number and type – longer incubations may be required for cells with weaker signal e.g. leukocytes.**
6. Measure the absorbance at 450 nm using a microplate reader. To measure the absorbance later, add 10 µl of 1% w/v SDS or 0.1 M HCl to each well, cover the plate and store it with protection from light at room temperature. No absorbance change should be observed for 24 hours.

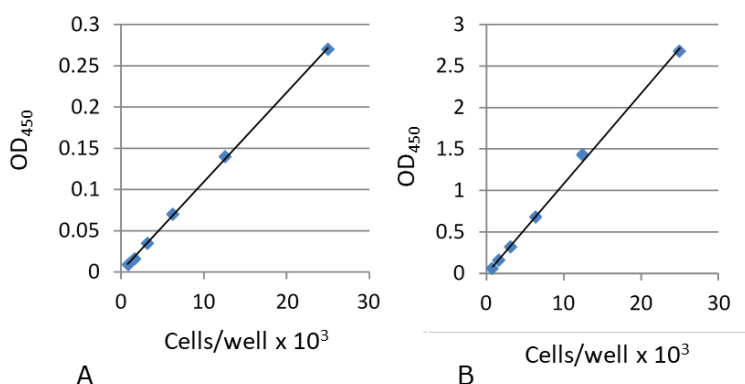


Figure 3. Standard curves for (A) HL60 and (B) HeLa cells.

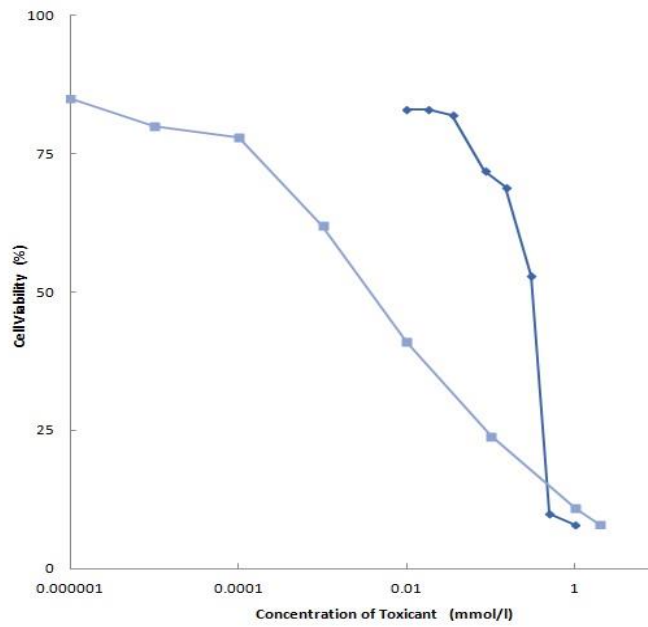


Figure 4. Toxicological test of chemicals using Orangu™.

Cell line:	HeLa
Culture medium:	MEM, 10% FBS
Chemicals:	Mitomycin-C (MMC) light blue Dodecylsulfate, sodium salt (SDS) dark blue
Incubation:	37°C, 5% CO <sub>2</sub> , 2 hours
Detection:	450 nm
Reference:	650 nm

## Notes on use of Orangu™

- The Orangu™ assay is based on the detection of dehydrogenase activity in viable cells. Conditions that affect dehydrogenase activity in viable cells may lead to a discrepancy between the actual viable cell number and the cell number determined using the Orangu™ assay.
- WST-8 may react with reducing agents to generate WST-8 formazan. Measure the background absorbance if reducing agents are used in cytotoxicity assays or cell proliferation assays.
- Take care not to introduce any bubbles into the wells – they may interfere with the absorbance reading.
- Culture media containing phenol red can be used with this kit for cell viability assays.
- Membrane filtration is recommended for the sterilization of the Orangu™ solution, if necessary.
- The incubation time varies by the type and number of cells in a well. For example, please see data in Figure 3 comparing HL60 and HeLa cells. Leukocytes give weak coloration, thus a long incubation time (up to 4 hours) or a large number of cells ( $\sim 1 \times 10^5$  cells/well) may be necessary.
- The cytotoxicity of this kit is very low. Further colour development is possible after reading the absorbance.
- Other cell proliferation assays, such as neutral red or crystal violet can be used after the Orangu™ assay on the same samples.
- Measure and subtract the absorbance at 600 nm or higher from that for a highly turbid cell suspension.



## Troubleshooting

### How many cells should there be in a well?

- For adherent cells, at least  $1 \times 10^3$  cells are necessary per well (100  $\mu$ l medium). For leukocytes, at least  $2.5 \times 10^3$  cells are necessary per well (100  $\mu$ l medium) because of low sensitivity. The recommended maximum number of cells per well for the 96-well plate is  $2.5 \times 10^4$ . If a 24-well or 6-well plate is used for this assay, calculate the number of cells per well accordingly, and adjust the volume of the Orangu™ solution in a well to 10% of the total volume.

### Is Orangu™ toxic to cells?

- Since the toxicity of Orangu™ is very low, the same cells can be used for other cell proliferation assays such as DNA fluorometric assays after the Orangu™ assay is completed.

### Does phenol red affect the assay?

- No. The absorption value of phenol red in a culture medium can be removed by subtracting the absorption value of a blank solution from the absorption value of each well. Therefore, a medium containing phenol red is usable for the Orangu™ assay.

### Does Orangu™ stain viable cells?

- No. Since WST-8 and its formazan dye are highly water-soluble, Orangu™ cannot be utilized for cell staining purpose.

### I do not have a 450 nm filter. What other filters can I use?

- You can use filters with an absorbance between 430 and 490 nm, even though 450 nm filter gives the best sensitivity.

## References

- M. Ishiyama et al. Talanta 44: 1299–1305. 1997
- M. Tominaga et al. Analytical Communications 36: 47-50. 1999



## Purchaser Notification

Limited warranty Cell Guidance Systems and/or its affiliate(s) warrant their products as set forth in the Terms of Sale found on the Cell Guidance Systems web site at [www.cellgs.com/Pages/Terms\\_and\\_Conditions.html](http://www.cellgs.com/Pages/Terms_and_Conditions.html)

If you have any questions, please contact Cell Guidance Systems.

This product incorporates licensed technologies. The purchase of this product conveys to the purchaser the limited, nontransferable right to use the purchased amount of the product only to perform internal research for the sole benefit of the purchaser. No right to resell this product or any of its components is conveyed. This product is for internal research purposes only and is not for use in commercial services of any kind, including, without limitation, reporting the results of purchaser's activities for a fee or other form of consideration. For information on obtaining additional rights, please contact [info@cellgs.com](mailto:info@cellgs.com).

CELL GUIDANCE SYSTEMS AND/OR ITS AFFILIATE(S) DISCLAIM ALL WARRANTIES WITH RESPECT TO THIS DOCUMENT, EXPRESSED OR IMPLIED, INCLUDING BUT NOT LIMITED TO THOSE OF MERCHANTABILITY, FITNESS FOR A PARTICULAR PURPOSE, OR NON-INFRINGEMENT. TO THE EXTENT ALLOWED BY LAW, IN NO EVENT SHALL CELL GUIDANCE SYSTEMS AND/OR ITS AFFILIATE(S) BE LIABLE, WHETHER IN CONTRACT, TORT, WARRANTY, OR UNDER ANY STATUTE OR ON ANY OTHER BASIS FOR SPECIAL, INCIDENTAL, INDIRECT, PUNITIVE, MULTIPLE OR CONSEQUENTIAL DAMAGES IN CONNECTION WITH OR ARISING FROM THIS DOCUMENT, INCLUDING BUT NOT LIMITED TO THE USE THEREOF.

Cell Guidance Systems' reagents and services enable control, manipulation and monitoring of the cell, both *in vitro* and *in vivo*.



**Growth Factors**

- Recombinant
- Sustained Release

**Exosomes**

- Purification
- Detection
- Tracking
- NTA Service

**Small Molecules**

**Cell Counting Reagent**

**Matrix Proteins**

**Cell Culture Media**

- Pluripotent Stem Cells
- Photostable
- *In Vitro* Blastocyst Culture
- ETS-embryo Culture
- Custom Manufacturing Service

**Gene Knock-Up System**

**Cytogenetics Analysis**



General [info@cellgs.com](mailto:info@cellgs.com)  
Technical Enquiries [tech@cellgs.com](mailto:tech@cellgs.com)  
Quotes [quotes@cellgs.com](mailto:quotes@cellgs.com)  
Orders [order@cellgs.com](mailto:order@cellgs.com)

[www.cellgs.com](http://www.cellgs.com)

**EUROPE**

**Cell Guidance Systems Ltd**  
Maia Building  
Babraham Bioscience Campus  
Cambridge  
CB22 3AT  
United Kingdom  
T +44 (0) 1223 967316  
F +44 (0) 1223 750186

**USA**

**Cell Guidance Systems LLC**  
Helix Center  
1100 Corporate Square Drive  
St. Louis  
MO 63132  
USA  
T 760 450 4304  
F 314 485 5424