

---

Product Manual


# OxiSelect™ Intracellular ROS Assay Kit (Green Fluorescence)

## Catalog Number

STA-342	96 assays
STA-342-5	5 x 96 assays

**FOR RESEARCH USE ONLY**  
Not for use in diagnostic procedures

---

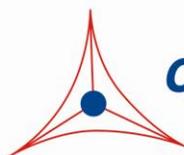

**CELL BIOLABS, INC.**  
*Creating Solutions for Life Science Research*

## **Introduction**

Accumulation of reactive oxygen species (ROS) coupled with an increase in oxidative stress has been implicated in the pathogenesis of several disease states. The role of oxidative stress in vascular diseases, diabetes, renal ischemia, atherosclerosis, pulmonary pathological states, inflammatory diseases, and cancer has been well established. Free radicals and other reactive species are constantly generated *in vivo* and cause oxidative damage to biomolecules, a process held in check by the existence of multiple antioxidant and repair systems as well as the replacement of damaged nucleic acids, proteins and lipids. Measuring the effect of antioxidant therapies and ROS activity intracellularly is crucial to suppressing or treating oxidative stress inducers.


Cell Biolabs' OxiSelect™ Intracellular ROS Assay Kit (Green Fluorescence) is a cell-based assay for measuring hydroxyl, peroxy, or other reactive oxygen species activity within a cell. The assay employs the cell-permeable fluorogenic probe 2', 7'-Dichlorodihydrofluorescein diacetate (DCFH-DA). In brief, DCFH-DA is diffused into cells and is deacetylated by cellular esterases to non-fluorescent 2', 7'-Dichlorodihydrofluorescein (DCFH), which is rapidly oxidized to highly fluorescent 2', 7'-Dichlorodihydrofluorescein (DCF) by ROS (Figure 1). The fluorescence intensity is proportional to the ROS levels within the cell cytosol. The effect of antioxidant or free radical compounds on DCFH-DA can be measured against the fluorescence of the provided DCF standard. The kit has a DCF detection sensitivity limit of 10 pM. Each kit provides sufficient reagents to perform up to 96 assays, including standard curve and unknown samples.

## **Assay Principle**

The OxiSelect™ Intracellular ROS Assay Kit is a cell-based assay for measuring antioxidant or ROS activity. Cells are cultured in a 96-well cell culture plate and then pre-incubated with DCFH-DA, which is cell-permeable (Figure 1). The unknown antioxidant or ROS samples are then added to the cells. After a brief incubation, the cells can be read on a standard fluorescence plate reader. The ROS or antioxidant content in unknown samples is determined by comparison with the predetermined DCF standard curve.

## **Related Products**

1. STA-320: OxiSelect™ Oxidative DNA Damage ELISA Kit (8-OHdG Quantitation)
2. STA-330: OxiSelect™ TBARS Assay Kit (MDA Quantitation)
3. STA-341: OxiSelect™ Catalase Activity Assay Kit
4. STA-343: OxiSelect™ Hydrogen Peroxide Assay Kit
5. STA-344: OxiSelect™ Hydrogen Peroxide/Peroxidase Assay Kit
6. STA-345: OxiSelect™ ORAC Activity Assay Kit
7. STA-347: OxiSelect™ In Vitro ROS/RNS Assay Kit (Green Fluorescence)
8. STA-350: OxiSelect™ Comet Assay Kit (3-Well Slides), 15 assays
9. STA-832: OxiSelect™ MDA Adduct Competitive ELISA Kit
10. STA-838: OxiSelect™ HNE Adduct Competitive ELISA Kit

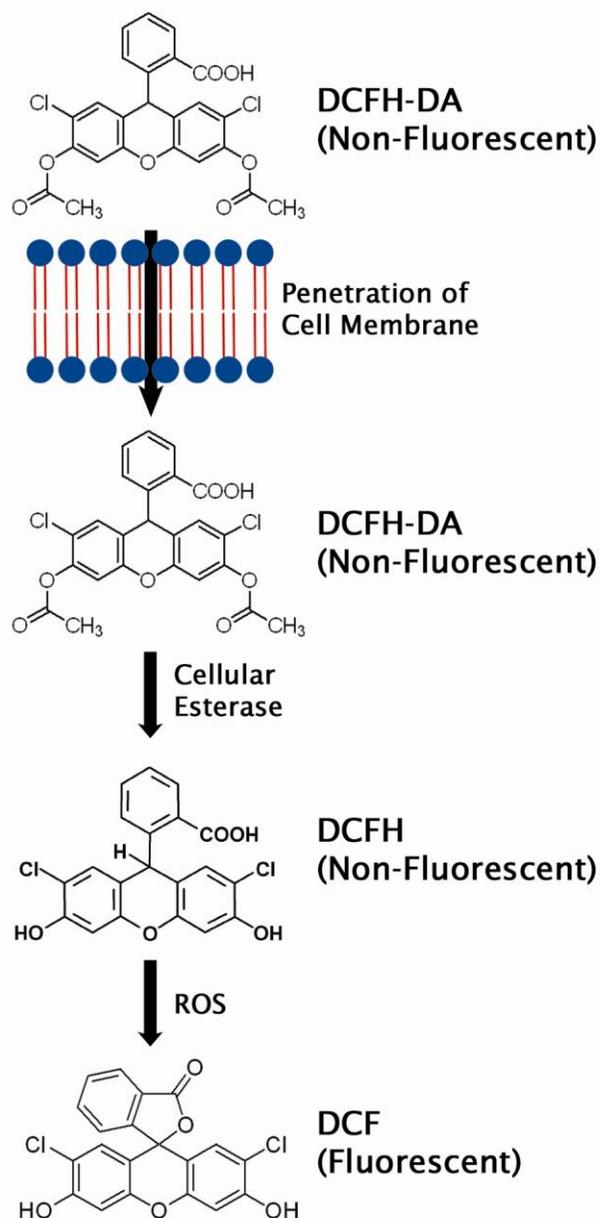

**Figure 1.** Mechanism of DCF Assay

### **Kit Components**

1. 20X DCFH-DA (Part No. 234201): One 500  $\mu$ L amber tube of a 20 mM solution in methanol.
2. DCF Standard (Part No. 234202): One 100  $\mu$ L amber tube of a 1 mM solution in DMSO.
3. Hydrogen Peroxide (Part No. 234102): One 100  $\mu$ L amber tube of an 8.821 M solution.
4. 2X Cell Lysis Buffer (Part No. 234203): One 20 mL bottle.

## **Materials Not Supplied**

1. Sterile DPBS for washes and buffer dilutions
2. Hank's Balanced Salt Solution (HBSS)
3. Cell culture medium (ie: DMEM +/-10% FBS)
4. 96-well black or fluorescence microtiter plate
5. Fluorescent microplate reader capable of reading 480 nm (excitation) and 530 nm (emission)

## **Storage**

Upon receipt, store the DCFH-DA and DCF Standard at -20°C. Avoid multiple freeze/thaw cycles. Store the Cell Lysis Buffer and Hydrogen Peroxide at 4°C.

## **Preparation of Reagents**

- 1X DCFH-DA: Dilute the 20X DCFH-DA stock solution to 1X in cell culture media, preferably without FBS. Stir or vortex to homogeneity. Prepare only enough for immediate applications.

*Notes:*

- *1X DCFH-DA/media solution contains 5% methanol. For cells that are sensitive to methanol, we recommend instead preparing a 0.1X (100 µM) solution of DCFH-DA in cell culture media.*
- *Due to light-induced auto-oxidation, DCFH-DA solutions at any concentration must be protected from light.*
- Hydrogen Peroxide (H<sub>2</sub>O<sub>2</sub>): Prepare H<sub>2</sub>O<sub>2</sub> dilutions in DMEM or DPBS as necessary. Do not store diluted solutions. Hydrogen Peroxide may be used as a positive control in the assay, or as a cell treatment.

## **Preparation of Standard Curve**

1. Prepare a 1:10 dilution series of DCF standards in the concentration range of 0 µM – 10 µM by diluting the 1 mM DCF stock in cell culture media (see Table 1).

<b>Standard Tubes</b>	<b>DCF Standard (µL)</b>	<b>Culture Medium (µL)</b>	<b>DCF (nM)</b>
1	10	990	10,000
2	100 of Tube #1	900	1000
3	100 of Tube #2	900	100
4	100 of Tube #3	900	10
5	100 of Tube #4	900	1
6	100 of Tube #5	900	0.1
7	100 of Tube #6	900	0.01
8	0	1000	0

**Table 1. Preparation of DCF Standards**

2. Transfer 75  $\mu\text{L}$  of each DCF standard to a 96-well plate suitable for fluorescence measurement. Add 75  $\mu\text{L}$  of the 2X Cell Lysis Buffer.
3. Read the fluorescence with a fluorescence plate reader at 480 nm excitation /530 nm emission.

## **Assay Protocol**

### **I. DCF Dye Loading**


1. Prepare and mix all reagents thoroughly before use. Each unknown sample should be assayed in duplicate or triplicate.
2. Culture cells in either a clear or black 96-well cell culture plate.  
*Note: If using a black plate, choose an appropriate plate based on your fluorometer's reader (i.e. choose a clear bottom black plate for bottom readers).*
3. Remove media from all wells and discard. Wash cells gently with DPBS or HBSS 2-3 times. Remove the last wash and discard.
4. Add 100  $\mu\text{L}$  of 1X DCFH-DA/media solution to cells. Incubate at 37°C for 30-60 minutes.
5. Remove solution. Repeat step three using multiple washes with DPBS or HBSS. Remove the last wash and discard.
6. Treat DCFH-DA loaded cells with desired oxidant or antioxidant in 100  $\mu\text{L}$  medium.

### **II. Quantitation of Fluorescence**


- Fluorescence microscopy or Flow cytometry: Fluorescence can be analyzed on an inverted fluorescence microscope or by flow cytometry using excitation and emission wavelengths of 480 nm and 530 nm, respectively.
- Fluorescence Plate Reader:
  - Assays performed in black cell culture fluorometric plates: Plate may be read immediately for kinetic analysis or after 1 hour for static analysis. Plates read for kinetic analysis may be read in increments of 1 and 5 minutes up to 1 hour or more as necessary. Read the fluorescence with a fluorometric plate reader at 480 nm/530 nm.
  - Assays performed in clear cell culture plates: After treatment with desired oxidant or antioxidant, carefully remove treatment media from all wells and discard. Wash cells gently with DPBS or HBSS 2-3 times. Remove the last wash and discard. Add 100  $\mu\text{L}$  of medium to each well. Add 100  $\mu\text{L}$  of the 2X Cell Lysis Buffer, mix thoroughly and incubate 5 minutes. Transfer 150  $\mu\text{L}$  of the mixture to a 96-well plate suitable for fluorescence measurement. Read the fluorescence with a fluorometric plate reader at 480 nm/530 nm.

## **Example of Results**


The following figures demonstrate typical ROS Assay results. Fluorescence measurement was performed on SpectraMax Gemini XS Fluorometer (Molecular Devices) with a 485/538 nm filter set and 530 nm cutoff. One should use the data below for reference only. This data should not be used to interpret actual results.

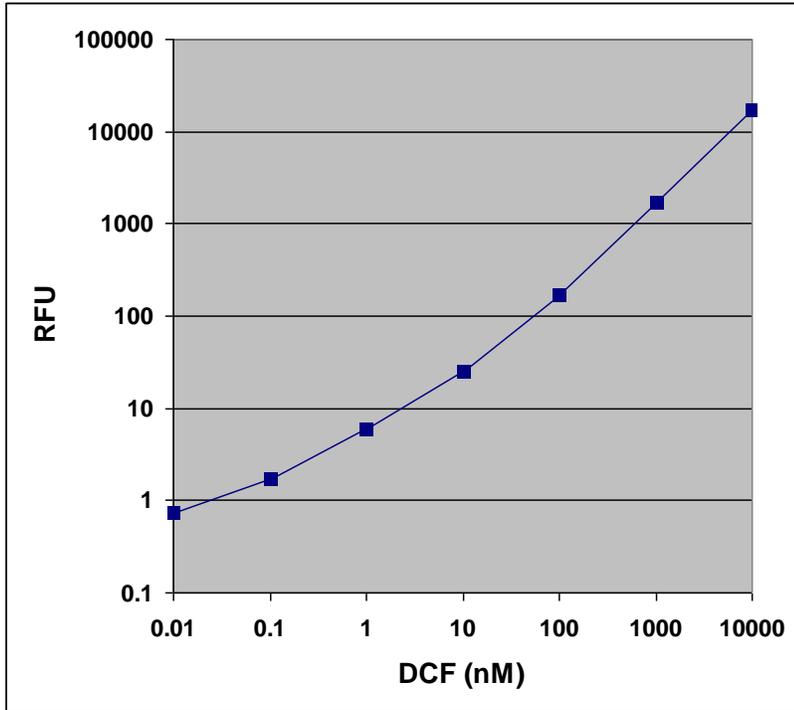

**Figure 2. DCF Standard Curve.**

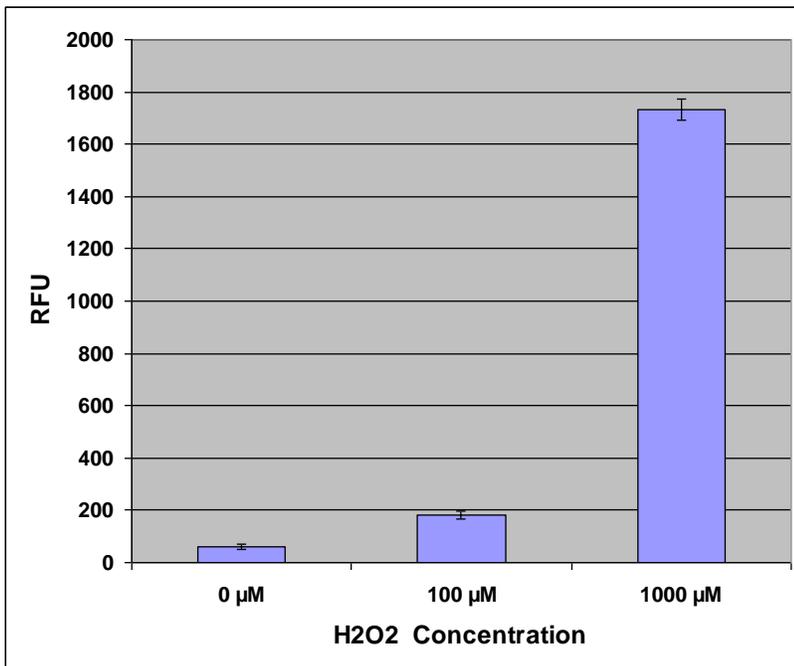

**Figure 3. ROS in HeLa cells treated with H<sub>2</sub>O<sub>2</sub>.** 50,000 HeLa cells in a 96-well plate were first pretreated with 1 mM DCFH-DA for 60 minutes at 37°C. Cells were then treated with various concentrations of H<sub>2</sub>O<sub>2</sub> for 20 minutes.

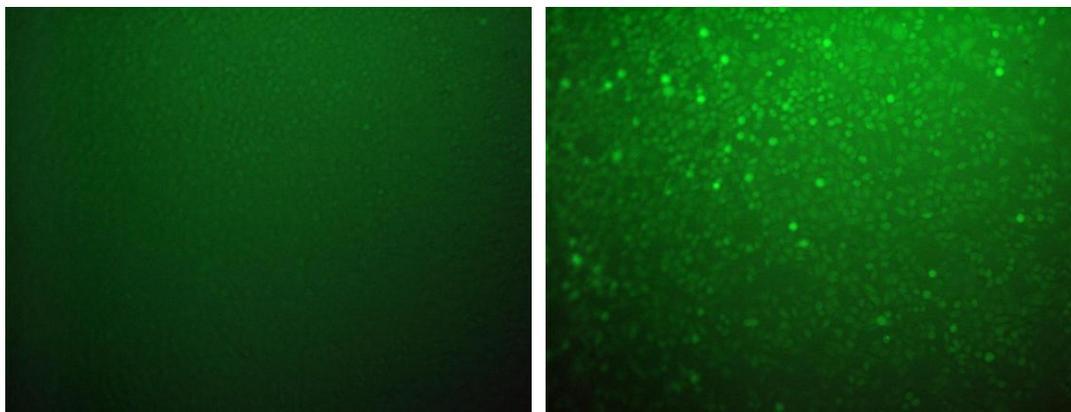

**Figure 4. DCF Fluorescence in H<sub>2</sub>O<sub>2</sub> treated HeLa cells after 1 hour. Left: 0 μM; Right: 1000 μM.**

### **References**

1. Bass DA, Parce JW, Dechatelet LR, Szejda P, Seeds MC, Thomas M. Flow cytometric studies of oxidative product formation by neutrophils: A graded response to membrane stimulation. *J Immunol.* 1983; **130**:1910-1917.
2. Brandt R, Keston AS. Synthesis of diacetyldichlorofluorescein: A stable reagent for fluorometric analysis. *Anal Biochem.* 1965; **11**:6-9.
3. Keston AS, Brandt R. The fluorometric analysis of ultramicro quantities of hydrogen peroxide. *Anal Biochem.* 1965; **11**:1-5.

### **Recent Product Citations**

1. Lall, R. K. et al. (2016). Dietary flavonoid fisetin increases abundance of high-molecular-mass hyaluronan conferring resistance to prostate oncogenesis. *Carcinogenesis*. doi:10.1093/carcin/bgw071.
2. Kobayashi, K. et al. (2016). Involvement of PARK2-mediated mitophagy in idiopathic pulmonary fibrosis pathogenesis. *J Immunol.* doi:10.4049/jimmunol.1600265.
3. Pan, Z. et al. (2016). β-Ecdysterone protects SH-SY5Y cells against 6-hydroxydopamine-induced apoptosis via mitochondria-dependent mechanism: Involvement of p38MAPK–p53 signaling pathway. *Neurotox Res.* doi:10.1007/s12640-016-9631-7.
4. Zhang, J. et al. (2016). Effect of macrophages on in vitro corrosion behavior of magnesium alloy. *J Biomed Mater Res Part A.* doi:10.1002/jbm.a.35788.
5. Ni, C. et al. (2016). Anesthetic isoflurane induces DNA damage through oxidative stress and p53 pathway. *Mol Neurobiol.* doi:10.1007/s12035-016-9937-8.
6. Crookenden, M. A. et al. (2016). Parturition in dairy cows temporarily alters the expression of genes in circulating neutrophils. *J Dairy Sci.* doi:10.3168/jds.2015-10877.
7. Behrens, J. et al. (2016). Coronin 1C-free primary mouse fibroblasts exhibit robust rearrangements in the orientation of actin filaments, microtubules and intermediate filaments. *Eur J Cell Biol.* doi:10.1016/j.ejcb.2016.04.004.
8. Chai, H. J. et al. (2016). Peptides from hydrolysate of lantern fish (*Benthosema pterotum*) proved neuroprotective in vitro and in vivo. *J Funct Foods.* **24**:438-449.
9. Tang, X. et al. (2016). Cystine deprivation triggers programmed necrosis in VHL-deficient renal cell carcinomas. *Cancer Res.* **76**:1892-1903.

10. Zhelev, Z. et al. (2016). Docosahexaenoic acid sensitizes leukemia lymphocytes to barasertib and everolimus by ROS-dependent mechanism without affecting the level of ROS and viability of normal lymphocytes. *Anticancer Res.* **36**:1673-1682.
11. Modepalli, N. et al. (2016). Transdermal delivery of iron using soluble microneedles: dermal kinetics and safety. *J Pharm Sci.* **105**:1196-1200.
12. Song, J. J. et al. (2016). Effect of cadmium on human middle ear epithelial cells. *J Int Adv Otol.* **11**:183-7.
13. Douard, N. et al. (2016). Impact of the chemical composition of poly-substituted hydroxyapatite particles on the in vitro pro-inflammatory response of macrophages. *Biomed Microdevices.* **18**:1-9.
14. Paul, A. et al. (2016). Nanoengineered biomimetic hydrogels for guiding human stem cell osteogenesis in three dimensional microenvironments. *J Mater Chem B.* doi:10.1039/C5TB02745D.
15. Gavalda-Navarro, A. et al. (2016). Changes in the expression of the human adenine nucleotide translocase isoforms condition cellular metabolic/proliferative status. *Open Biol.* doi:10.1098/rsob.150108.
16. Subramani, R. et al. (2016). Nimbolide inhibits pancreatic cancer growth and metastasis through ROS-mediated apoptosis and inhibition of epithelial-to-mesenchymal transition. *Sci Rep.* doi:10.1038/srep19819.
17. Tu, G. et al. (2016). Allicin attenuates H<sub>2</sub>O<sub>2</sub>-induced cytotoxicity in retinal pigmented epithelial cells by regulating the levels of reactive oxygen species. *Mol Med Rep.* doi:10.3892/mmr.2016.4797.
18. Johnson, R. et al. (2016). Aspalathin, a dihydrochalcone C-glucoside, protects H9c2 cardiomyocytes against high glucose-induced shifts in substrate preference and apoptosis. *Mol Nutr Food Res.* doi:10.1002/mnfr.201500656.
19. Yamaguchi, Y. et al. (2016). Arsenic acid inhibits proliferation of skin fibroblasts, and increases cellular senescence through ROS mediated MST1-FOXO signaling pathway. *J Toxicol Sci.* **41**:105-113.
20. Kuranaga, Y. et al. (2016). Anti-oncogenic gem-dihydroperoxides induce apoptosis in cancer cells by trapping reactive oxygen species. *Int J Mol Sci.* doi:10.3390/ijms17010071.
21. Liu, Y. et al. (2016). Mefloquine effectively targets gastric cancer cells through phosphatase-dependent inhibition of PI3K/Akt/mTOR signaling pathway. *Biochem Biophys Res Commun.* doi:10.1016/j.bbrc.2016.01.046.
22. Maronedze, C. et al. (2016). Cyclic nucleotide monophosphates in plants and plant signaling. doi:10.1007/164\_2015\_35.
23. Li, C. et al. (2015). The interplay between autophagy and apoptosis induced by tanshinone IIA in prostate cancer cells. *Tumor Biol.* doi:10.1007/s13277-015-4602-9.
24. Di, K. et al. (2015). Marizomib activity as a single agent in malignant gliomas: ability to cross the blood-brain barrier. *Neuro Oncol.* doi:10.1093/neuonc/nov299.
25. Slusarczyk, J. et al. (2015). Anti-inflammatory properties of tianeptine on lipopolysaccharide-induced changes in microglial cells involve Toll-like receptor-related pathways. *J Neurochem.* doi:10.1111/jnc.13452.
26. Zhelev, Z. et al. (2015). 2-deoxy-D-glucose sensitizes cancer cells to barasertib and everolimus by ros-independent mechanism (s). *Anticancer Res.* **35**:6623-6632.
27. Sayes, C. M. et al. (2015). Physicochemical characteristics of two prototypical home-use consumer products containing engineered nanomaterials. *J Environ Anal Toxicol.* **5**:2161-0525.

28. Yu, C. L. et al. (2015). Ferulic acid protects against lead acetate-induced inhibition of neurite outgrowth by upregulating HO-1 in PC12 cells: Involvement of ERK1/2-Nrf2 pathway. *Mol Neurobiol.* doi:10.1007/s12035-015-9555-x.
29. Gao, S. et al. (2015). Apoptotic effects of Photofrin-Diomed 630-PDT on SHEEC human esophageal squamous cancer cells. *Int J Clin Exp Med.* **8**:15098.
30. Chung, E. H. et al. (2015). miR-505 downregulates 6-phosphofructo-2-kinase/fructose-2, 6-biphosphatase 4 to promote cell death in glioblastoma. *Cancer Transl Med.* **1**:158-165.
31. Wang, Y. et al. (2015). ROS-mediated activation of JNK/p38 contributes partially to the pro-apoptotic effect of ajoene on cells of lung adenocarcinoma. *Tumor Biol.* doi:10.1007/s13277-015-4181-9.
32. Kim, K. A. & Yim, J. E. (2015). Antioxidative activity of onion peel extract in obese women: a randomized, double-blind, placebo controlled study. *J Cancer Prev.* **20**:202-207.
33. Kim, Y. et al. (2015). Insecticidal activity of chlorine dioxide gas by inducing an oxidative stress to the red flour beetle, *Tribolium castaneum*. *J Stored Prod Res.* **64**:88-96.
34. Ha, Y. R. et al. (2015). Effect of farnesyltransferase inhibitor R115777 on mitochondria of *plasmodium falciparum*. *Korean J Parasitol.* **53**:421-30.
35. Hikita, H. et al. (2015). Activation of the mitochondrial apoptotic pathway produces reactive oxygen species and oxidative damage in hepatocytes that contribute to liver tumorigenesis. *Cancer Prev Res (Phila).* **8**:693-701.
36. Dong, H. et al. (2015). Paeoniflorin inhibition of 6-hydroxydopamine-induced apoptosis in PC12 cells via suppressing reactive oxygen species-mediated PKC $\delta$ /NF- $\kappa$ B pathway. *Neuroscience.* **285**:70-80.

Please see the complete list of product citations: <http://www.cellbiolabs.com/reactive-oxygen-species-ros-assay>.

## **Warranty**

These products are warranted to perform as described in their labeling and in Cell Biolabs literature when used in accordance with their instructions. THERE ARE NO WARRANTIES THAT EXTEND BEYOND THIS EXPRESSED WARRANTY AND CELL BIOLABS DISCLAIMS ANY IMPLIED WARRANTY OF MERCHANTABILITY OR WARRANTY OF FITNESS FOR PARTICULAR PURPOSE. CELL BIOLABS' sole obligation and purchaser's exclusive remedy for breach of this warranty shall be, at the option of CELL BIOLABS, to repair or replace the products. In no event shall CELL BIOLABS be liable for any proximate, incidental or consequential damages in connection with the products.

## **Contact Information**

Cell Biolabs, Inc.  
 7758 Arjons Drive  
 San Diego, CA 92126  
 Worldwide: +1 858-271-6500  
 USA Toll-Free: 1-888-CBL-0505  
 E-mail: [tech@cellbiolabs.com](mailto:tech@cellbiolabs.com)  
[www.cellbiolabs.com](http://www.cellbiolabs.com)

©2009-2016: Cell Biolabs, Inc. - All rights reserved. No part of these works may be reproduced in any form without permissions in writing.