

How Did My Cells Die?

Choosing the Right Apoptosis Assay



Apoptosis encompasses a non-inflammatory cascade of molecular events and results in distinct morphological changes including reduction in cellular volume, nuclear fragmentation, condensation of chromatin, DNA degradation, cell membrane blebbing, and subsequent dispersal into apoptotic bodies. Due to the variety of apoptotic stimuli and subtleties in resulting pathways, it can be critical to determine which correlates of cell death are affected by your experimental conditions.

Mitochondrial Membrane Potential


During apoptosis, the potential of the proton gradient across the mitochondrial inner membrane decreases in association with the opening of the mitochondrial permeability pore and the release of apoptogenic factors, such as cytochrome *c*. In some apoptotic models, loss of membrane potential is thought to be an early event in the apoptotic process. However, others suggest that loss of membrane potential is a consequence of the apoptotic signaling pathway. A simple way to assess membrane potential of mitochondria in a cell population is the use of positively charged dyes, such as JC-1 and TMRE, that accumulate in the electronegative interior of active mitochondria.

JC-1/TMRE

JC-1 changes from red fluorescence when aggregated in healthy mitochondria to green fluorescence when membrane potential is lost, as in apoptosis. TMRE accumulates in polarized mitochondria and is suitable for monitoring changes in mitochondrial membrane potential in live cells. Depolarized mitochondria with decreased membrane potential are unable to accumulate TMRE.

Caspase Activation

Caspases are activated *via* an extrinsic pathway when extracellular death ligands (such as FasL or TNF- α) bind to transmembrane death receptors, recruiting initiator caspase-8. Caspases are also activated intrinsically by cytochrome *c* released from mitochondria. In response, an apoptosome assembles to recruit and

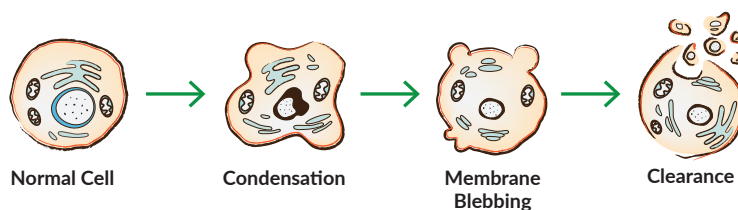

activate initiator caspase-9. Initiator caspases activate executioner caspases-3 and -7, which cleave specific substrates that directly result in the morphological changes that define apoptosis.

Caspase-3/7

The protease activity of caspase-3 and -7 can be detected by using the sequence-specific fluorogenic substrate N-Ac-DEVD-N'-MC-R110, which, upon cleavage, generates a fluorescent product.

Nuclear Fragmentation, Chromatin Condensation, and DNA Degradation

Nuclear fragmentation results from the disintegration of the nuclear lamina *via* proteolysis by caspases and the collapse of the nuclear envelope. Another distinguishing characteristic of apoptosis is the condensation of chromatin accompanied by the hydrolysis of nuclear DNA into a ladder of fragments. Hoechst 33342, a cell-permeable, fluorescent nucleic acid stain, is often used to identify nuclear condensation by microscopic analysis. The Golgi, endoplasmic reticulum, and mitochondrial networks also undergo fragmentation during apoptosis. As they are dismantled, the pieces are dispersed into plasma membrane blebs.

DNA Stains

Hoechst dyes are cell permeable and bind to nucleic acids in live or fixed cells. Their blue emission spectrum is convenient for researchers who plan to use multiple fluorescent stains in a single sample. An alternate cell-permeable DNA stain is DRAQ5™, whose far-red excitation can be multiplexed with fluorophores with blue to orange emission. Membrane-impermeant dyes, such as propidium iodide, DAPI, DRAQ7™, or RedDot™2, are ideal for specifically staining the nuclei of dead cells, whose plasma membrane integrity is compromised.

Membrane Blebbing

Because of cytoskeleton weakening, the cell cytoplasm pushes against unsupported areas of the plasma membrane. Membrane protrusions, termed blebs, can be visualized microscopically. This is thought to result from myosin-dependent contraction of cortical bundles of actin, pushing the cytosol against the cell cortex and causing blebs in areas where the cytoskeleton has deteriorated. Membrane blebbing is a key step toward apoptotic cells budding off into smaller apoptotic bodies.

Apoptotic Clearance Signals

Early in apoptosis, apoptotic cells secrete “find-me” signals to recruit engulfing cells and display “eat-me” signals that engulfing cells can recognize. This select interaction between apoptotic cells and engulfing cells ensures specific targeting of dying cells for non-inflammatory clearance. “Find-me” signals that recruit phagocytes include lysophosphatidylcholine, sphingosine-1-phosphate, fractalkine, and nucleotides like ATP and UTP. Through caspase-mediated channel opening, plasma membrane pannexin 1 (PANX1) channels release nucleotide “find-me” signals to recruit phagocytes. Caspase-dependent PANX1 channel opening also allows the entry of a small group of fluorescent monomeric cyanine dyes including TO-PRO®-3. Thus, pannexin channel functioning can be taken advantage of to identify cells that are sending “find-me” signals.

TO-PRO®-3

TO-PRO®-3 is a far-red fluorescent dye that enables detection of a very early event in the initiation of apoptosis. It enters cells through selective pannexin channels that are activated in a caspase-dependent manner early in the apoptotic process. Thus, cells that might be considered viable due to lack of detection of traditional apoptosis markers (*i.e.*, Annexin V) can be identified as having just begun the apoptotic process.

Once the phagocyte is attracted to the apoptotic cell, it recognizes phosphatidylserine and phosphatidylethanolamine residues redistributed from the inner to outer leaflet of the phospholipid bilayer that are now exposed on the cell surface. This membrane rearrangement—a well-recognized hallmark of apoptosis—designates the apoptotic cell

as different from its viable counterparts and, thus, is a sign for phagocytes to eat the foreign cell. The phospholipid binding protein Annexin V adheres to phosphatidylserine residues exposed during apoptosis. It is used to probe for phosphatidylserine on the outer membrane of apoptotic cells in order to identify the “eat-me” phase of apoptosis involving the redistribution of membrane phospholipids.

Annexin V

Annexin V can be conjugated to various fluorochromes (*e.g.*, FITC, PE, APC) to probe for phosphatidylserine on the outer membrane of apoptotic cells. Staining for Annexin V can be combined with markers of plasma membrane integrity (*i.e.*, RedDot™2, propidium iodide, DAPI) to distinguish apoptotic cells from necrotic cells.

Multiplex is Best

These distinguishing morphological features and key parts of the apoptotic signaling pathway can be observed and measured to verify the mode of cell death in your model system. However, a single readout can be misleading under some experimental conditions (*e.g.*, distinguishing apoptosis from necrosis). Multiplexed assays are available to provide a more complete understanding of the apoptotic process in a single sample. For example, measuring TO-PRO®-3, Annexin V, TMRE, and DAPI allows you to concomitantly evaluate “find-me” *versus* “eat me” signals, mitochondrial membrane potential, and nuclear fragmentation for a more complex picture of how apoptosis is affecting your cells. Early apoptosis markers (*i.e.*, TO-PRO®-3), enable cells that might have been considered viable due to a lack of traditional apoptosis markers (*i.e.*, Annexin V) to now be identified as having just begun the apoptotic process. Thus, a multiplexed method that includes multiple markers for various stages of apoptosis allows for the quantitative segmentation of the total sample population into distinct phases: viable, early apoptosis, late apoptosis, apoptotic bodies, and non-cellular debris.

The range of cell-based assays available from Cayman is designed to give you the flexibility and efficiency needed to determine the variable stages of cell death in your model system.

Kit Recommendations

Apoptosis Assays

Item No.	Product Name	Measure	Readout	Time to Results
601410	Annexin V APC Assay Kit	Phosphatidylserine residues on apoptotic cells ("eat-me" signal) Includes DAPI to assess plasma membrane permeability	Flow cytometer	30 minute staining protocol
600300	Annexin V FITC Assay Kit	Phosphatidylserine residues on apoptotic cells ("eat-me" signal) Includes propidium iodide to assess plasma membrane permeability	Flow cytometer Fluorescence microscope	30 minute staining protocol
601420	Annexin V PE Assay Kit	Phosphatidylserine residues on apoptotic cells ("eat-me" signal) Includes DAPI to assess plasma membrane permeability	Flow cytometer	30 minute staining protocol
10009135	Caspase-3/7 Fluorescence Assay Kit	Caspase-3/7 enzymatic activity levels in cell lysates	Fluorescence plate reader	2 hours
601360	Early Apoptosis Detection Assay Kit	TO-PRO®-3: Caspase-activated PANX1 ("find-me" signal) Annexin V FITC: Phosphatidylserine residues on apoptotic cells ("eat-me" signal) TMRE: Mitochondrial membrane potential DAPI: Plasma membrane permeability	Flow cytometer	30 minute staining protocol
10009172	JC-1 Mitochondrial Membrane Potential Assay Kit	Mitochondrial membrane potential	Flow cytometer Fluorescence microscope Fluorescence plate reader	<1 hour staining protocol
701560	JC-1 Mitochondrial Membrane Potential Flow Cytometry Assay Kit	Mitochondrial membrane potential	Flow cytometer	15 minute staining protocol
601430	Mitochondrial PT Pore Assay Kit	Calcein AM/Cobalt chloride: Inner mitochondrial membrane integrity TMRE: Mitochondrial membrane potential Includes controls for mitochondrial membrane potential depolarization (FCCP) and calcein quenching	Flow cytometer Fluorescence microscope	15 minute staining protocol
601280	Multi-Parameter Apoptosis Assay Kit	Annexin V FITC: Phosphatidylserine residues on apoptotic cells ("eat-me" signal) TMRE: Mitochondrial membrane potential RedDot™2: Plasma membrane integrity Hoechst 33342: Nuclear morphology	Flow cytometer Fluorescence microscope	<1 hour staining protocol
701310	TMRE Mitochondrial Membrane Potential Assay Kit	Mitochondrial membrane potential	Fluorescence plate reader	<1 hour staining protocol

Related Assays

Cell Proliferation Assays

Item No.	Product Name	Measure	Readout	Time to Results
10009349	Cell Cycle Phase Determination Kit	Cell cycle progression	Flow cytometer	~3 hours
10009853	CFSE Cell Division Assay Kit	Up to eight cell divisions in activated lymphocytes	Flow cytometer	<2 hour staining protocol
10009365	MTT Cell Proliferation Assay Kit	Cell viability and proliferation in cultured cells	Colorimetric plate reader	8-22 hours
10008883	WST-1 Cell Proliferation Assay Kit	Cell viability and proliferation in cultured cells	Colorimetric plate reader	2-4 hours
10010199	WST-8 Cell Proliferation Assay Kit	Cell viability and proliferation in cultured cells	Colorimetric plate reader	2-4 hours
10010200	XTT Cell Proliferation Assay Kit	Cell viability and proliferation in cultured cells	Colorimetric plate reader	~4 hours

Cytotoxicity Assays

Item No.	Product Name	Measure	Readout	Time to Results
600120	7-AAD/CFSE Cell-Mediated Cytotoxicity Assay Kit	Cytotoxicity in immune effector cells	Flow cytometer	6-7 hours
700410	ATP Detection Assay Kit – Luminescence	Total ATP levels in cell lysates	Luminescence plate reader	<30 minutes
600140	Autophagy/Cytotoxicity Dual Staining Kit	Regulation of autophagy and cytotoxicity in cultured cells	Fluorescence microscope Fluorescence plate reader	20 minute staining protocol
601170	LDH Cytotoxicity Assay Kit	Drug-induced cytotoxicity in cultured cells	Colorimetric plate reader	<2 hours

Metabolism Assays

Item No.	Product Name	Measure	Readout	Time to Results
700410	ATP Detection Assay Kit - Luminescence	Total ATP levels in cell lysates	Luminescence plate reader	<30 minutes
701220	Calcium Assay Kit	Calcium in plasma, serum, urine, tissue homogenates, and cell lysates	Colorimetric plate reader	<30 minutes
10007640	Cholesterol Fluorometric Assay Kit	Total cholesterol in plasma or serum	Fluorescence plate reader	<1 hour
10009582	Glucose Colorimetric Assay Kit	Glucose in plasma, serum, and urine	Colorimetric plate reader	<1 hour
700480	Glycogen Assay Kit	Glycogen from tissue samples	Fluorescence plate reader	<1 hour
700510	L-Lactate Assay Kit	L-Lactate in serum, plasma, whole blood, urine, saliva, and cell culture	Fluorescence plate reader	<1 hour
601470	LDL Uptake Flow Cytometry Assay Kit	LDL cellular uptake	Flow cytometer	~24 hours
700470	Pyruvate Assay Kit	Pyruvate in serum, plasma, blood, urine, and saliva	Fluorescence plate reader	<1 hour
10010303	Triglyceride Colorimetric Assay Kit	Triglyceride levels in plasma, serum, cell lysates, and tissue homogenates	Colorimetric plate reader	<1 hour
700320	Uric Acid Assay Kit	Uric acid in serum, plasma, and urine	Fluorescence plate reader	<1 hour

Standardization & Monitoring Assays

Item No.	Product Name	Measure	Readout	Time to Results
700460	Creatinine (serum) Colorimetric Assay Kit	Creatinine levels in plasma and serum	Colorimetric plate reader	<1 hour
500701	Creatinine (urinary) Colorimetric Assay Kit	Creatinine levels in urine	Colorimetric plate reader	<1 hour
10012445	His-Tag Detection ELISA Kit	His-tagged proteins from cell lysates or affinity column fractions	Colorimetric plate reader	<3 hours
10009325	Malachite Green Phosphate Assay Kit	Inorganic free phosphate in aqueous solutions	Colorimetric plate reader	<1 hour
701410	Protease Activity Assay Kit	Protease activity in samples	Fluorescence polarization plate reader Fluorescence plate reader	<30 minutes
704002	Protein Determination Kit	Total protein concentration	Colorimetric plate reader	<1 hour

Have a question for our Technical Support team?

Our technical support and product development scientists are here to help you choose the most appropriate assay for your experimental endpoint, so you can obtain the most accurate answers for your research questions.

Call us at [\(734\) 975-3888](tel:(734)975-3888)

Email us at techserv@caymanchem.com

Visit us online at www.caymanchem.com/TechnicalServices

View a complete list of our assay kits at www.caymanchem.com